

J.K.

## UAB "LT Biotech"

302303586, PVM kodas LT 100004741118,  
Rugijų 21-24, LT-08419, Vilnius Reg. Nr. 127918, V.Į.  
Registru centras Vilniaus filialas  
Tel/fax +370 5 216 02 27

### Fetal Bovine Serum, IgG Stripped

#### Product Description

Serum is the blood component that can be obtained after coagulation and by removing cellular components. Besides serum proteins it contains, e.g. growth factors, amino acids and hormones. This comprehensive mix makes serum one of the most important supplements, supporting cells to grow and proliferate *in vitro* cell culture.


Of special interest is the fetal bovine serum (FBS). It is especially rich in growth factors and is particularly low in antibodies, which may influence the cell culture work.

FBS is processed through protein A-Sepharose columns to reduce the IgG content of serum to below 5 µg/ml. This process is gentle to the product and ensures an unimpaired growth promotion. The specially treated FBS has a very low IgG antibody concentration and is outstandingly suitable for hybridoma cell culture. Thereby, the IgG stripped FBS facilitates the purification of monoclonal antibodies.

Product	Origin	Volume	Cat. No.
Fetal Bovine Serum IgG Stripped	Collected in South America	100 ml	LTfBS-SA100IGD
		500 ml	LTfBS-SA500IGD

#### Product Specifications

pH	6.8 – 8.2
Osmolality	240 – 340 mOsm/kg
Endotoxin	As reported
Total Protein	3.0 – 4.5 g/dl
IgG	< 5 µg/ml
Albumin	As reported
Hemoglobin	As reported
Mycoplasma	Not detected
Virus tested for	PI-3, BVDV, BVDV-AB, BHV-1
Sterility	Tested
Storage	Store at ≤-15°C

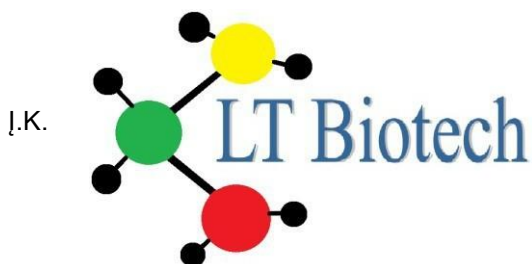

J.K.

## UAB “LT Biotech”

302303586, PVM kodas LT 100004741118,  
Rugių 21-24, LT-08419, Vilnius Reg. Nr. 127918, V.Į.  
Registrijų centras Vilniaus filialas  
Tel/fax +370 5 216 02 27

### **Additional optional treatments**

#### ***Heat inactivation***

Heat inactivation will inactivate the complement system, antibodies and other active enzymes. It has to be done in a carefully controlled process in order to avoid damaging the cell growth promoting properties of the serum and reducing the formation of unwanted precipitates.

The process involves heating the serum in a shaking water bath at exactly +56°C for 30 minutes. The shaking will help avoid the formation of protein and other forms of precipitates. After 30 minutes the serum is then cooled back down to room temperature as quickly as possible to avoid excessive exposure to heat which can damage e.g. growth factors and vitamins.

#### ***Gamma irradiation***

Capricorn Scientific has established process parameters and controls for maximum inactivation of contaminants by an innovative gamma irradiation process in small sized boxes at 25 – 35kGy. When FBS bottles are arranged in a pallet during gamma irradiation, high irradiation doses (58 kGy) are necessary to irradiate the centrally located bottles with the required dose (>30 kGy). This and the associated higher temperatures may affect the serum quality. If the irradiation is performed in single boxes, the maximum dose of irradiation is performed in single boxes, the maximum dose of irradiation does not exceed 38.7 kGy. Gentle irradiation as used by Capricorn Scientific is less likely to affect final serum quality.

#### **Quality control**

Only sera batches which pass our strict quality control are released for sale. Standard parameters which are determined include pH, osmolality, content of protein, albumin, IgG and hemoglobin, endotoxin level, sterility, mycoplasma detection and virus testing.

#### **Precautions and Disclaimer**

This product is for research use only. Please consult the Material Safety Data Sheet for information regarding hazards and safe handling practices.