

Lepta MMA HR

1. Basic product information

Lepta MMA HR is a modified agarose-based multimodal chromatography resin. Different from traditional ion exchange chromatography resin, it is composed of multiple groups with different properties: amino groups, hydrogen bonds and hydrophobic groups. The chromatographic resin can remove major impurities such as nucleic acid, host protein, aggregates and viruses, and is widely used in the fields of polysaccharides, antibodies, viral vectors or vaccines. Compared with traditional single-mode chromatography resin, Lepta MMA HR shows better performance:

- (1) It has high dynamic capacity, combining with most impurities in one step, saving time and cost.
- (2) The improved Lepta base frame is more rigid, so it can achieve higher process flow rate at lower back pressure and improve process efficiency.
- (3) It can tolerate a wide range of pH and conductivity, which is helpful to explore the optimal process conditions.
- (4) It has a fine particle size design and high resolution.

2. Chromatography resin parameters


Resin type	Multimodal
Functional group	Amino groups, hydrogen bonds and hydrophobic groups
Matrix	Highly cross-linked agarose
Median particle size	40 µm
Ion load	0.08–0.11 mmol Cl ⁻ /ml
Recommended flow rate	90–500 cm/h
Maximum flow rate	600 cm/h
Withstand pressure	5 bar
Use temperature	4–30°C

3. Chemical resistance

pH stability*	2–14
Chemical stability	Common aqueous solution, 30% isopropanol**, 75% ethanol**, 1M NaOH, 1M acetic acid, 6M guanidine hydrochloride, 8M urea
Avoid	Oxidising agents, anionic detergents

* The physical and chemical properties and functions of the chromatographic resin have no obvious change after being placed in an environment of 40°C and pH 2–14 for 7 days.

** v/v, volume ratio

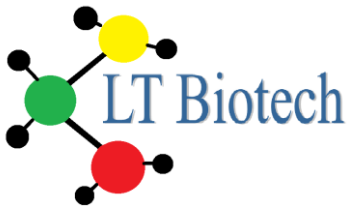

4. Method of use

4.1 Chromatographic conditions

(1) Buffer selection: a buffer salt whose buffer group does not interact with the chromatographic resin should be selected. If the binding and elution mode is used, the equilibration buffer should be a buffer with a high pH (usually 1 pH unit higher than the isoelectric point of the target molecule) to facilitate the binding of the target molecule, and the stability of the sample in the buffer should also be considered; the elution buffer is usually a buffer in which a high concentration of salt (such as 1M NaCl) is added to the equilibration buffer. If the flow-through mode is used, the equilibration buffer should adopt conditions that are conducive to the binding of impurities. After the target molecule has completely flowed through, it should be washed directly with high-concentration salt.

(2) Flow rate: generally choose a linear flow rate of 90–500 cm/h according to the height of the column bed.

(3) Sample pretreatment: In order to prevent the sample from clogging the column, the sample needs to be filtered with a 0.45 µm microporous membrane before loading the sample. It is recommended that the pH and conductivity of the sample are adjusted to be consistent with the equilibrium buffer (dilution, ultrafiltration can be used and desalting to adjust the pH and conductivity of the sample).

4.2 Chromatography steps

(1) Equilibration: use equilibration buffer to fully equilibrate the chromatography column until the pH and conductivity are stable and basically consistent with the equilibration buffer. This step usually requires 3–5 times the column volume (CV).

(2) Sample loading*: according to the binding capacity measured in the small test, determine the sample loading volume and loading amount of the sample on Lepta MMA HR.

(3) Impurity washing*: use equilibration buffer or other suitable buffer to wash the chromatography column until the UV stabilises and returns to the baseline.

(4) Elution*: elution is achieved by increasing the concentration of salt ions. The concentration of salt ions in the eluent can be gradually increased through a linear gradient or a step gradient to elute molecules with different binding strengths. pH gradient elution or mixed elution can also be used.

(5) Regeneration: rinse the column with a high-salt buffer (such as 2M NaCl).


(6) Re-equilibration: re-equilibrate the column with equilibration buffer.

*Note: If the flow-through mode is used, the sample loading step should be set to collect; the washing step, to ensure that the target molecules have completely flowed through, then stop the collection; the elution step should directly use high-salt buffer, and impurities can be washed away.

5. Cleaning and regeneration

Contaminants (e.g. lipids, endotoxins and proteins) accumulate on the column as the number of uses of the chromatography resin increases. Determine the frequency of CIP according to the degree of contamination of the chromatography resin (if the contamination is considerable, CIP is recommended after each use to ensure repeatability of results and to prolong the working life of the chromatography resin). For different types of impurities and contaminants, the recommended cleaning conditions are as follows:

- Removal of strongly binding proteins: wash with 5 CV of 2M NaCl solution or use a high salt buffer not lower than pH 2, such as 1M NaAc solution.
- Removal of strongly hydrophobic proteins and precipitated proteins: first wash with 5 CV of 1M NaOH solution, then wash the lye with 5–10 CV of ultra pure or pure water.

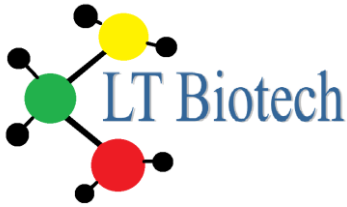

- Removal of lipoproteins and lipids: first wash with 5 CV of 70% ethanol or 30% isopropanol, then rinse with 5–10 CV of ultra pure or pure water.

Note: 70% ethanol or 30% isopropanol should be degassed before use; the flow rate should be 30–60 cm/h during CIP. Reverse cleaning should be used when the clogging is severe.

To reduce the microbial load, it is recommended that 0.5~1M NaOH solution is used to treat the chromatography resin for 15~30 minutes.

6. Storage

Keep the unopened chromatography resin in the original container and store at 4~30°C in a well-ventilated, dry and clean place. Do not freeze. Wash the used column with 2–3 CV of 20% ethanol solution and store at 2~8°C.

7. Destruction and recycling

Since chromatography resin is difficult to degrade in nature, it is recommended that the waste chromatography resin is incinerated to protect the environment. For chromatography resin that has been in contact with biologically active samples such as viruses and blood, follow the local biosafety requirements before destroying or disposing of it.

8. Packing method

Detailed information on resin packaging is available on request. Please contact your local distributor.

9. Ordering information

Product name: Lepta MMA HR

Product Cat. No	Package
613-00025	25 ml
613-00100	100 ml
613-00500	500 ml
613-01000	1 L
613-05000	5 L
613-10000	10 L
613-20000	20 L